

STANDARD OPERATING PROCEDURES

DIVISION OF COMPARATIVE MEDICINE
UNIVERSITY OF SOUTH FLORIDA

SOP#: 1161.2

Date Issued: 5/15

Date Revised: 7/20

Page 1 of 2

TITLE: Germinator™ 500 Glass Bead Dry Sterilizer
SCOPE: Animal Care Personnel
RESPONSIBILITY: Facility Manager, Professional and Administrative Staff
PURPOSE: To Outline the Proper Procedures for Use and Maintenance of a Germinator™ Glass Bead Sterilizer

I. PURPOSE

1. This procedure outlines the use and maintenance of the Germinator™ Glass Bead Sterilizer used to

